



Leuther laboratories

N 1006 County Road M ~ Coon Valley, WI 54623 ~ 608-788-8180

THIO-BAG®

Instructions for Whirl-Pak® THIO-BAGS®

Caution: Do not allow the thio tablet to fall out of the bag.

OPENING



Tear off top along perforation.



Pull tabs outward to open bag. Sometimes a pull on the bottom is also helpful.



Fill bag to required fill line. The 4 oz. fill line will ensure that 100 ml are collected.

CLOSING



Pull wire ends to close bag.



Whirl bag three complete revolutions. (Do not roll tapes down.) Leave airspace for mixing in lab.



Turn tape inward on opposite face of fold.

NASCO • FORT ATKINSON, WISCONSIN • MODESTO, CALIFORNIA • SPECTRUM/NASCO • NEWMARKET, ONTARIO, CANADA
PRINTED IN U.S.A. U05352 /RV 9-09

1. Remove screens, hoses and/or any other attachments on sample faucet.
2. Clean faucet inside and outside with regular bleach.
3. Rinse and run water for 5 minutes so no bleach remain in sample. Any sanitizers in sample will invalidate test.
4. Take sample using the Thio-Bag® as above, or sterile bottle provided by laboratory.
5. Keep sample at 40° F or less. Bring to lab on ice.
6. Sample must be tested with 30 hours to be a legal test.
7. Come by laboratory or contact us and we can mail Thio-Bags to you with instructions.